

# ReadySeq

## Preparation of samples



Sample is a sequencing reaction prepared by using the BigDye Terminator v3.1 Cycle Sequencing kit (Applied Biosystems), cleaned and finally dissolved in at least 15  $\mu$ l of Hi-Di formamide (Applied Biosystems). Samples must be delivered in 96-well plates (Standard type, not Fast), compatible with Applied Biosystems sequencers.

We recommend that immediately after finishing your sample preparation you send them on ice by using a courier service (mail or our drop boxes are not recommended, ready sequencing reactions must be stored on ice and analyzed fast).

**Samples must be delivered in 96-well plates. Do not use parafilm.**

We recommend you contact us before sending your first batch of samples!

v20210511

# ReadySeq